TRUWEIGH-2

Instruction Manual

Thank you for purchasing TW2 SERIES pocket scale.

Model: TW2-100 & TW2-600

Please read all operating instructions carefully before use. This electronic scale is a precision instrument. With normal care and proper treatment, it will provide years of reliable service.

Manual- English

NOTE:

- *Never load the scale with more than the maximal capacity.
- *Avoiding any exposure to extreme heat or cold, your scale works better when operated at normal room temperature. Always allow the scale to acclimate to normal room temperature for at least one hour before use.
- *Give your scale sufficient warm up time. Usually 30 to 60 seconds before calibration to give the internal components a chance to stabilize.
- *Keep your scale in a clean environment. Dust, dirt, moisture, vibration, air currents and a close distance to other electronic equipment can all cause an adverse effect on the reliability and accuracy of your scale.
- *Handle with care. Gently apply all items to be weighed onto tray top.
- Although this scale is designed to be quite durable, try to avoid rough treatment as this may permanently damage the internal sensor.
- *Avoid shaking, dropping or otherwise shocking the scale. This is a precision instrument and must be handled with extreme care.
- *Only operate the scale on a stable, vibration free surface.

OPERATION

Place scale on a horizontal flat surface

1. Press the [ON/T] key to turn the scale on.
2. Wait until "0.0" is displayed.
3. Place the object(s) on the weighing platform.
4. By pressing the [M] key, you can switch between weighing units.
5. To turn the scale off, press and hold the [ON/T] key.

II. TARE WEIGHING

1. Turn on the scale as described above.
2. Place a tray or container on the scale platform.
3. Press [ON/T] key and wait until "0.0" is displayed
4. Add the "net-weight-item".

III. CALIBRATION

1. Turn on the scale.
 2. Press and hold the [M] key until the LCD display "CAL", release the [M] key.
 3. Press [M] key again, the display will show "CAL" followed by the required calibration weight value.
 4. place the required calibration weight on the center of the platform, the display will show "PASS" then return to the normal weighing mode.
- Calibration is complete! You can turn off the scale.

IV SPECIFICATIONS

Capacity: 100g/1500gn/500ct/64.24dwt
Readability: 0.01g/0.15gn/0.05ct/0.01dwt
Units: g/gn/ct/dwt

Capacity: 600g/21.16oz/19.29ozt/385.8dwt
Readability: 0.1g/0.01oz/0.01ozt/0.1dwt
Units: g/oz/ozt/dwt

Tare Range: Tare full capacity
Auto Off: 180 seconds off
Operation temp: 10 to 30 degree C
Display: LCD 5 1/2 digits
Power: 2 x AAA batteries

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